

Biological Microscope

Model Number:STM-2030B/T

Introduction Manual



This instruction manual is for STM-2030B Binocular and STM-2030T Trinocular biological microscope. To insure safety and obtain optimum performance and familiarize yourself fully with the use of this microscope. We recommend that you read the manual thoroughly before operating the microscope, Attain this manual instruction in an easily accessible place near the microscope for the further reference.

CONTENTS

| | |
|---------------------------------|----|
| User Notice..... | 2 |
| 1.Components..... | 4 |
| 2.Assembly..... | 5 |
| 2.1 Assembly Diagram | 5 |
| 2.2 Assembly Steps | 6 |
| 3. Adjustment & Operation | 10 |
| 3.1 Adjustment Sets..... | 10 |
| 3.2 Operation..... | 11 |
| 4. Specification Table | 14 |
| 5. Outfit | 15 |
| 6. Troubleshooting Guide | 16 |

User Notice

I . Safety Notes


1. Carefully open the box, avoid the accessories, like lens, dropping to ground and being damaged.
2. Do keep the instrument out of direct sunlight, high temperature or humidity, dusty and easy shaking environment. Make sure the stage is smooth, horizontal and firm enough.
3. When moving the instrument, please use two hands to grip with the two sides of the microscope body.
4. When running, the lamp house and nearby parts will be very hot. Please ensure there is enough cooling room for them.
5. Make sure the instrument is earthed, to avoid lighting strike.
6. For safety, be sure the main switch is in “O”(off) state before replacing the halogen lamp or the fuse, then cut off the power, and do the operation after the lamp bulb and the lamp house completely cool down.(**Specified: Halogen Lamp6V/20W**)
7. Check the input voltage: be sure the input voltage which signed in the back of the microscope is consistent with the power supply voltage, or it will bring a serious damage to the instrument.
8. Use the factory supplied power cord, please.

II. Maintenance

1. All the lenses have been well checked and adjusted. It is forbidden to disassemble them yourself.
2. The nosepiece and coarse/fine focus unit have a compact and precise frame; please don't disassemble them as possibly as you can.
3. The nosepiece and coarse/fine focus unit have a compact and precise frame; please don't disassemble them as possibly as you can.
4. Keep the instrument clean, wipe dust regularly, and be attention to avoid contaminating the optical elements especially.
5. The contaminations on the prism, as finger mark and oil, could be gently wiped with a piece of soft cloth or tissue paper, gauze which has been immersed in pure alcohol or aether. (**Note that the alcohol and ether are highly flammable, do keep them away from the fire or potential sources of electrical sparks, and use them in a drafty room as possible as you can.**)
6. Do not attempt to use organic solvents to clean the microscope components other than the glass components. To clean them, use a lint-free, soft cloth slightly moistened with a diluted neutral detergent.

7. When using, if the microscope is splashed by liquid, cut off the power at once, and wipe up the moisture.
8. Do not disassemble any parts of the microscope, which will affect the function or decline the performance of the microscope.
9. Place the instrument in a cool, dry position. When not using the microscope, keep it covered with a dust cover. Make sure the lamp socket is cool before covering the microscope.

1.Components

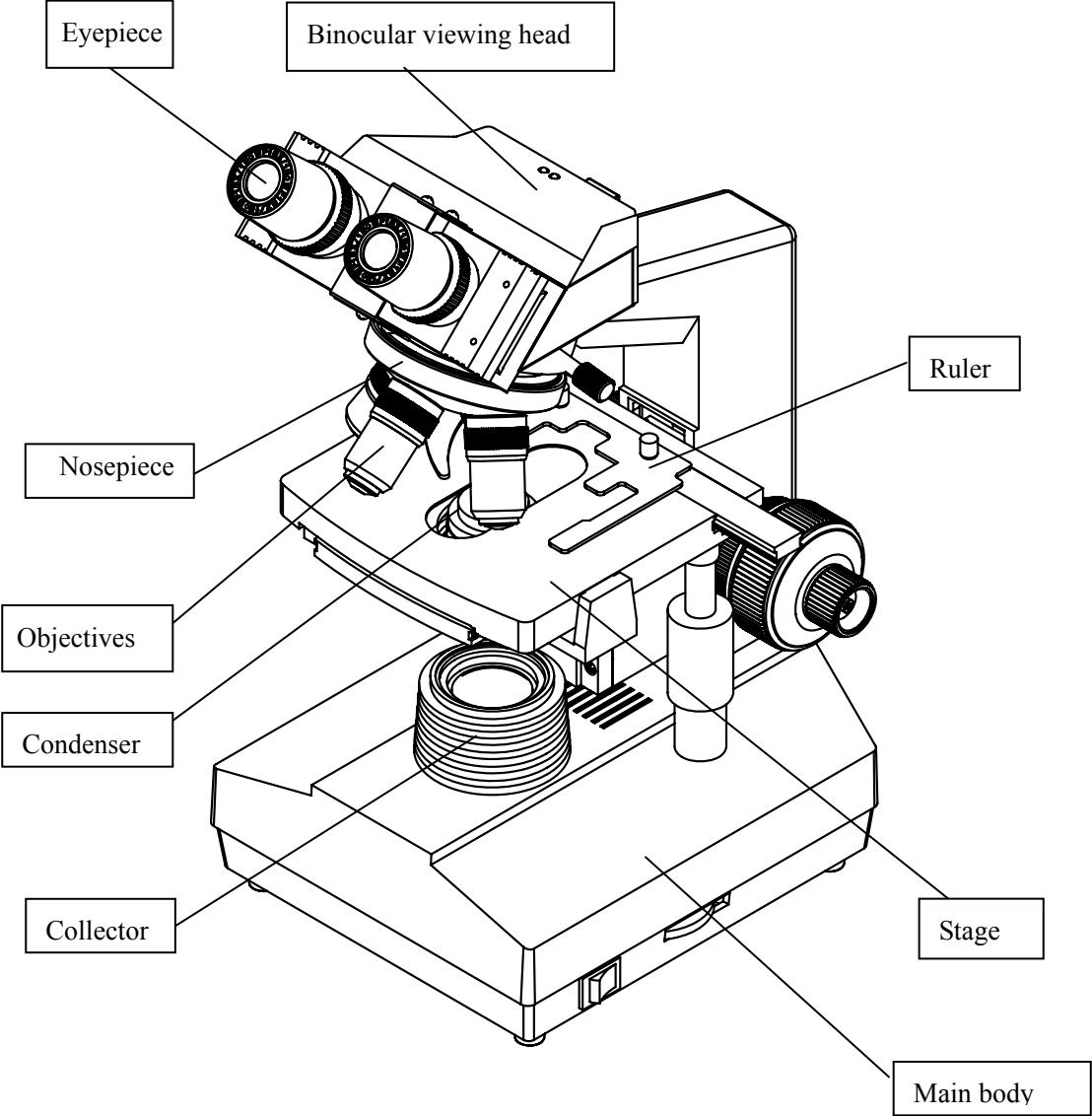


STM-2030B Biological Microscope

2.Assembly

2.1 Assembly Diagram

The following figure shows the installation sequence of the components. The number in the figure shows the assembly steps.

- ★ Before installing, be sure every components is clean, do not score any parts or glass surface.
- ★ Keep well with hexagon wrench provided. When replacing the components, you will need it again.

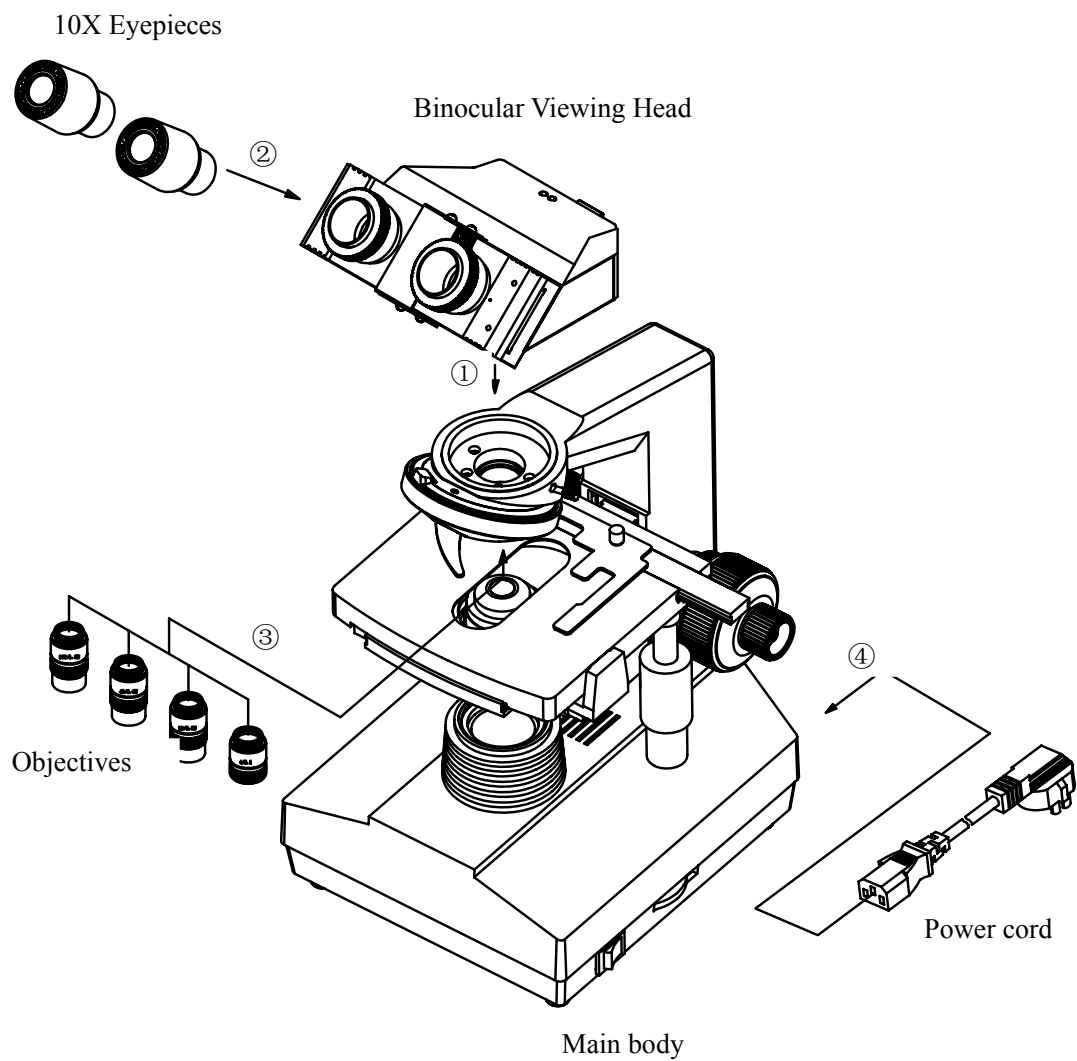

2.2 Assembly Steps

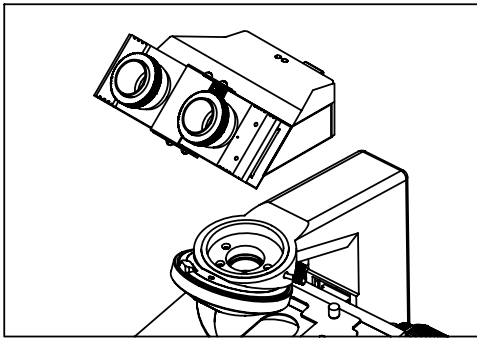


Fig.1

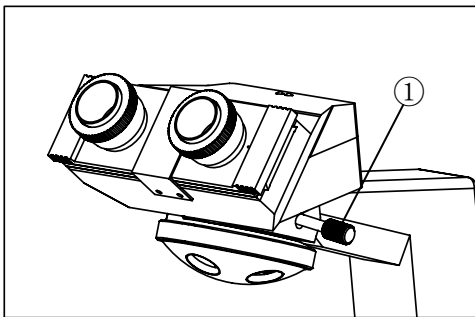


Fig.2

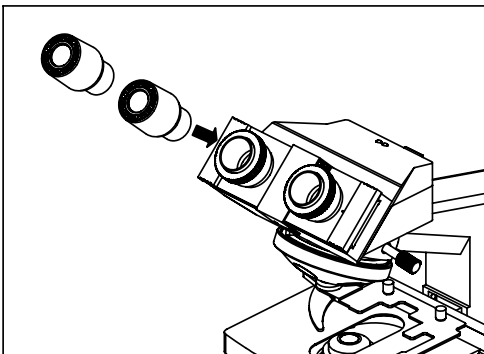


Fig.3

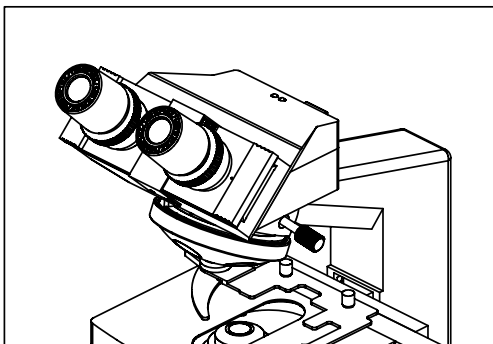


Fig.4

2.2.1 Installing Binocular viewing head (Fig.1, 2)

Insert the digital viewing head into the microscope head, turn into the right position, then screw down the bolt① to fix it.

2.2.2 Installing the eyepieces (Fig.3, Fig.4)

Insert the eyepieces into the eyepiece tube until they are against each other as shown in Fig.4.

NOTE:

Operation Conditions:

1. Temperature: $0^{\circ}\text{C} \sim 40^{\circ}\text{C}$, Maximum Relative Humidity: 85%.
2. High Temperature: High Temperature and humidity will result in a mildewing, dew and even ruinous instrument.
3. Avoid placing the instrument in a dusty environment. When ending your microscope operation, please cover it with the dust cap.
4. Lay the microscope in a plan and stable position, please.

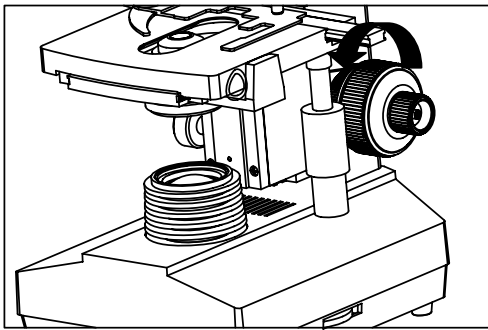


Fig.5

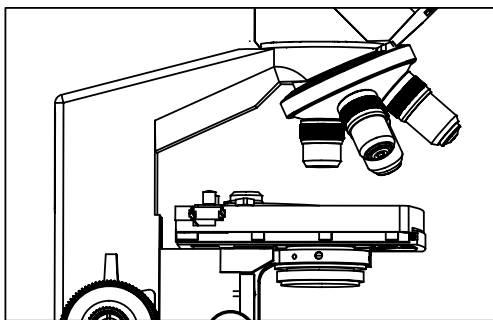


Fig.6

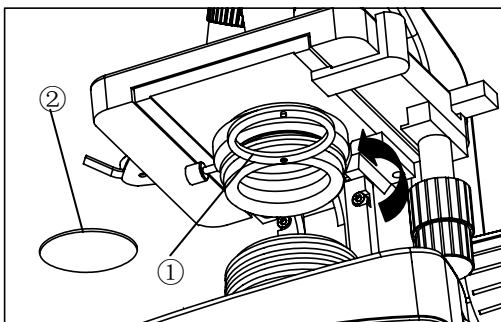


Fig.7

2.2.3 Installing objectives (Fig.5& 6)

1. Adjusting the coarse focus knob until the support device of the mechanical stage reaches its low limit position.
2. Screw the lowest magnification objective into the nosepiece from the left or the right side, then revolve the nosepiece clockwise and mount other objectives by the sequence of low to high magnification

◇ Installing objective this way will make the change of magnification to be easier during using.

- ★ Clean the objective regularly, for lens is susceptible to dust.
- ★ When operating, use 10×magnification objective to search and focus specimen firstly, then replace with higher magnification objective if necessary.
- ★ When replacing the objective, slowly turn the nosepiece until you hear “clicked”, which means the objective is in the required position--the light path center.

2.2.4 Installing the color filters (Fig.7)

1. Turn the condenser bracket① out at the direction of arrow in Fig.7
2. Put the required filters② into the holder on the bracket, and then turn the bracket back to the right position.

★ Baby blue and green filters are available in standard outfit.

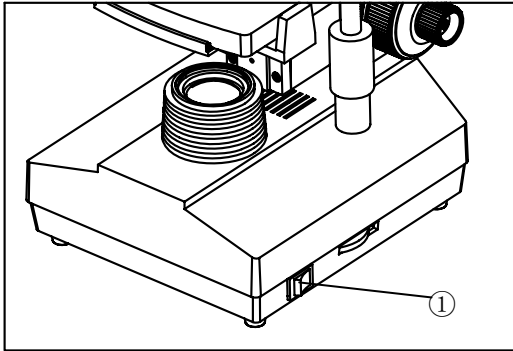


Fig.8

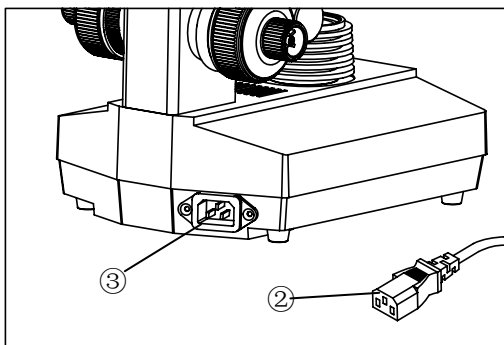


Fig.9

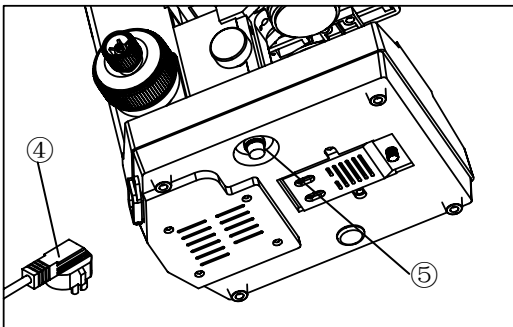


Fig.10

2.2.5 Connecting the power cord (Fig.8, 9,10)

★ The cable and cords are vulnerable when bent or twisted, never subject the power cord to excessive force.

1. Turn the main switch① to “O” (off) state before connecting the power cord.
2. Insert the power plugs② into the power jack③ of the microscope; make sure the connection is well.
3. Plug the power cord ④ into the power supply receptacle safely. Make sure the connection is well.

★ Do use the supplied power cord all the time. If lost or damaged, select the same standard cord, please.

2.2.6 Replacing the Fuse (Fig.8, 9, 10)

Do remember to set the main switch① to the state of “O” (OFF) and unplug the power cord ②before replacing the fuse. Rotate the fuse kits⑤ out of the holder, replace with a new fuse, then rotate it back to the holder again.

★ The rating of the new fuse should be the same as the old one.

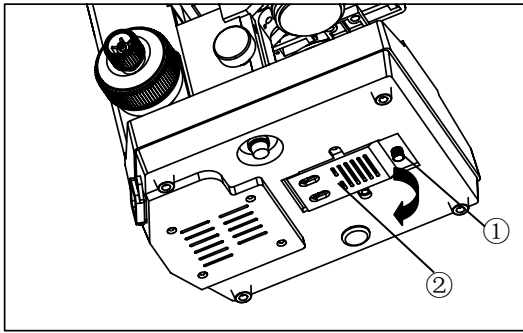


Fig.11

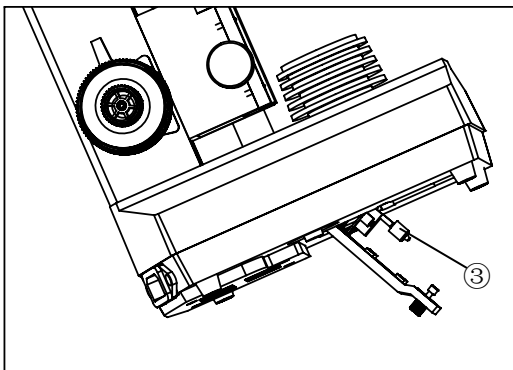


Fig.12

2.2.7 Installing and replacing the lamp (Fig.11, 12)

✧ Please use the specified halogen Lamp 6V20W. Replace the lamp as follows if necessary.

1. Please set the main switch to “O” (off) state before replacing, and make sure the bulb, the lamp room and periphery are all cool enough to carry no burn. Then, you can do your replacing.
2. Loose the bolt①and open the window② on the bottom of the microscope base with “—”type screwdriver.
3. Pull out the old bulb③, hold the new bulb after you wrap it with gauze or other protection materials and insert its pin as deeply as possible into the jack in the lamp holder.
4. Close the window and tighten the bolt①.

★ Please insert the bulb gently, or it will be damaged by excessive extrusion.

★ Do not touch the halogen bulb with bare hands. It will shorten the service life or cause it to burst. If you leave fingerprints on the surface carelessly, clean it with a piece of dry soft cloth.

3. Adjustment & Operation

3.1 Adjustment Sets (Fig.14, Fig.15)

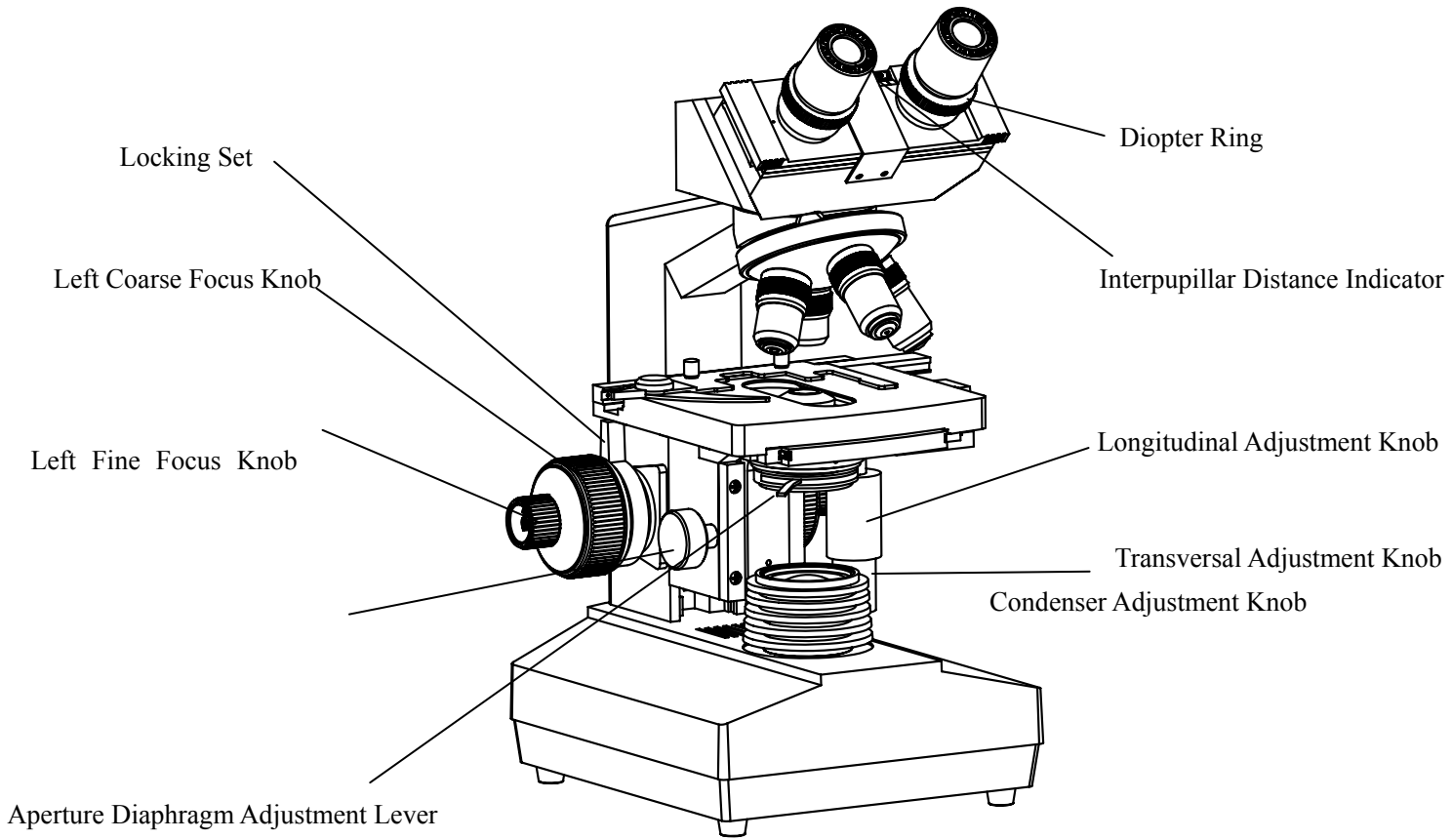


Fig.14

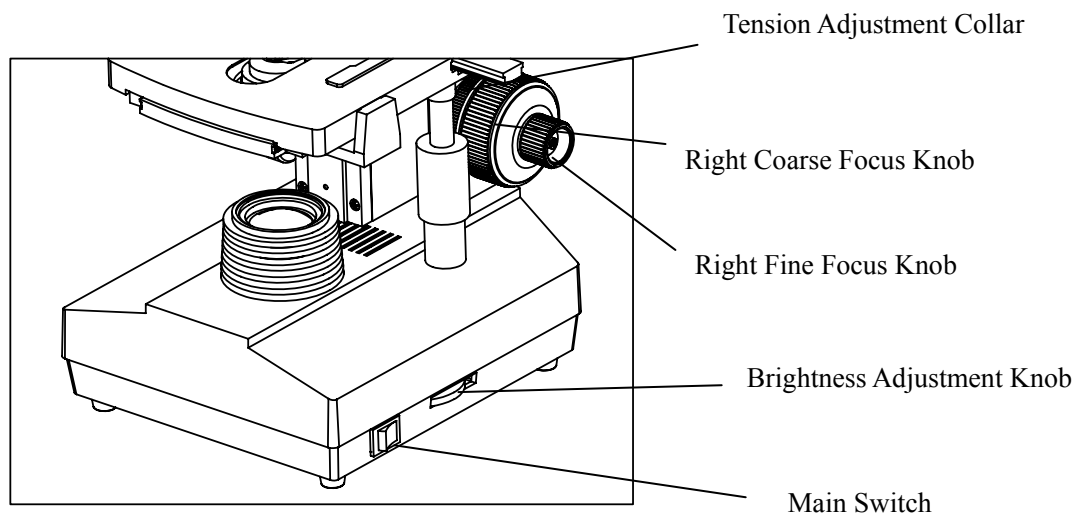


Fig.15

3.2 Operation

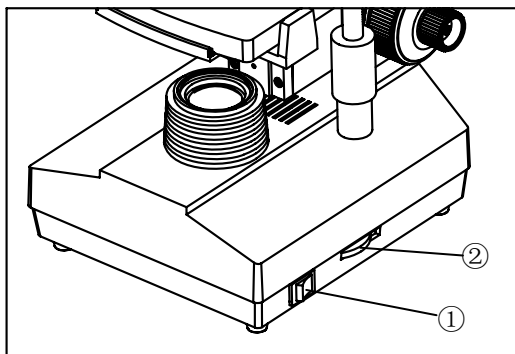


Fig.16

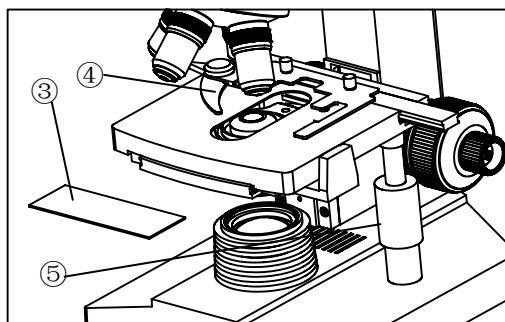


Fig.17

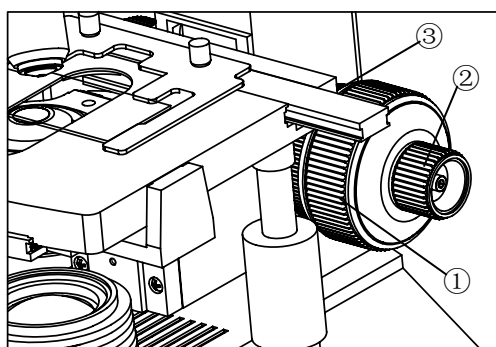


Fig.18



Fig.19

3.2.1 Adjusting the brightness (Fig.16)

1. Connect the power, turn on the main switch①(shown in the figure)which on the bottom side of the base to “—”(on).
2. Turning the brightness adjustment knob ②clockwise, the voltage decline, and the brightness weaken; Whereas turning at the opposite direction, the voltage raise, and the brightness strengthen.

★ Using the microscope at a lower voltage can prolong the service life of the bulb.

3.2.2 Placing the specimen (Fig.17)

1. Place the specimen③ on the center of the stage, and then nip it with the specimen holder④.
2. Turn the transversal and longitudinal adjustment knobs which on the mechanical ruler to move the specimen onto the required position.

★ Be careful when changing the objective. If you finish the observation with the short working distance objective, and want to change another one, be careful of not letting the objective touch the specimen.

3.3.3 Focusing the specimen (Fig.18, 19)

1. Focus the specimen with 10X objective. To avoid the objective touching the specimen during focusing, you should raise the mechanical stage to let the specimen close to the objective at first, then slowly separate them to bring the specimen to focus.
2. Turn the coarse focus knob ①conversely to lower the specimen and search images in the 10×ocular simultaneously, and then use the fine knob② to make focus. After that, you can replace with other magnification objectives safely, and focus without the risk of damaging the specimen.
3. When the specimen is focused, tighten the locking set (in Fig.14) to give a height limit to the stage. In this way, the stage only can lower, avoiding false operation and damage to the specimen.

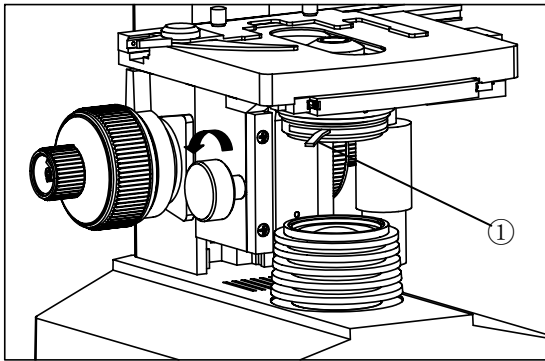


Fig.20

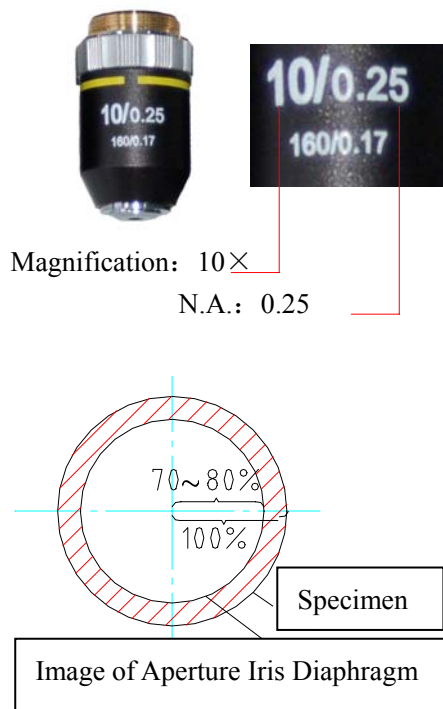


Fig.21

- The tight tension of the coarse focus knob has already been adjusted before leaving factory. If loosen (e.g. the stage slip down by its weight), please screw the intention adjustment collar③ to the right position by the supplied spanner.

3.3.4 Condenser Adjustment (Fig.20)

Turn the condenser focus knob to move the condenser up and down. Raise the condenser when using the high magnification objective, and descend it when using the low magnification one.

- ★ The condenser and the objective are coaxial. It has been adjusted before leaving factory, so the user needn't to adjust them by self (the distance between the top of the condenser and the stage should be in the range of 0.03mm~0.4mm.)
- ★ The highest position of the condenser has been adjusted too. It also needn't any user's operation.

3.3.5 Aperture Iris Diaphragm Adjustment (Fig.20,21)

Turn the aperture iris diaphragm lever① to adjust the aperture iris diaphragm.

Generally, setting the aperture iris diaphragm to 70- 80% of the N.A. of the objective in use will provide an image with good contrast.

- If the size of the aperture diaphragm minified, the brightness and the resolution declined, while the contrast and the depth of field increased; In other words, if the size largen, the brightness and the resolution improved, but the contrast and the depth of field declined.
- Generally, setting the size of the condenser aperture diaphragm at 70%~80% of the numerical aperture, you can obtain a clear image with enough contrast. If the open of the aperture diaphragm is too small, the resolution were very low, so please don't

- Minify the aperture below 60% of the objective's numerical aperture unless in a special case, for instance, observing an almost transparent specimen.
- The numerical aperture is marked on the objective. For example, the mark “ 10/0.25 ” means the magnification is 10×, and the numerical aperture is 0.25.
- If you want to observe the image of the aperture iris diaphragm, remove one eyepiece and look through the tube. You will see a dark circle encroaching on the bottom of the tube.

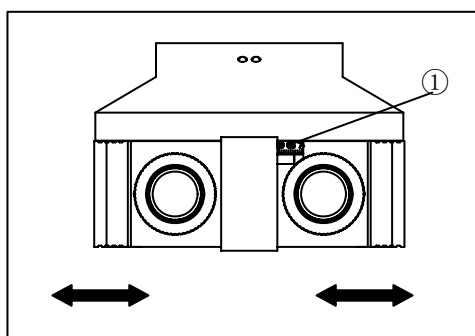


Fig.22

3.3.6 Adjusting the Interpupillary Distance (Fig.22)

The interpupillary distance range:

55mm~75mm.

While looking through the eyepieces, move both sliders to adjust interpupillary distance until the left and right fields of view coincide completely. The value on the right slider is your interpupillary distance.

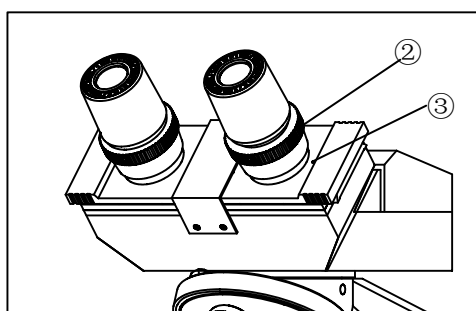


Fig.23

3.3.7 Adjusting the Diopter (Fig.23)

Turn the diopter rings② on both eyepieces to align the number on the rings (which is corresponding to your interpupillary distance) with the white dot③ on both sliders to bring the specimen into focus.

★ The diopter range of the eyepiece is ± 5 diopter. The number aligned to the line on the viewing head is the diopter in use.

4. Specification Table

4.1 Main specifications

| | |
|------------------------|---|
| Mechanical Tube Length | 160mm |
| Viewing Head | Sliding Binocular/Trinocular Head, 45°Inclined, Interpupillary Distance 55-75 mm |
| Eyepiece | Field of view: ϕ 18mm |
| Nosepiece | External Quadruple Nosepiece |
| Objective | Achromatic objectives 4 \times , 10 \times , 40 \times , 100 \times (immersion oil) |
| Focusing | Coaxial Coarse & Fine Adjustment, Moving Range 30mm, Fine Division 0.004mm. |
| Condenser | Abbe Condenser, NA=1.25 |
| Stage | Double Layers Mechanical Stage 140mm \times 140mm, Moving Range 74 \times 50mm |
| Illumination | LED illumination or 6V/20W Halogen Lamp |

4.2 Eyepieces and Objectives

1. Achromatic Objectives

| Magnification | Numerical Aperture (NA) | Thickness of glass slide (mm) | Focal length (mm) | Working Distance(mm) | Type |
|---------------|-------------------------|-------------------------------|-------------------|----------------------|------|
| 4 \times | 0.10 | 0.17 | 31.05 | 18 | Dry |
| 10 \times | 0.25 | 0.17 | 17.13 | 6.5 | Dry |
| 40 \times | 0.65 | 0.17 | 4.65 | 0.53 | Dry |
| 100 \times | 1.25 | 0.17 | 2.906 | 0.13 | Oil |

Plan Achromatic Objectives are optional

| Designation | Magnification | Numerical Aperture | Working Distance |
|---------------------------|---------------|--------------------|------------------|
| Plan Achromatic Objective | 4X | 0.10 | 15.8mm |
| | 10X | 0.25 | 12.2 mm |
| | 40X(S) | 0.65 | 0.37 mm |
| | 100X(Oil,S) | 1.25 | 0.13 mm |

1. Eyepieces

| Designation | Magnification | Field of view(mm) | Focal Length(mm) |
|-------------|---------------|-------------------|------------------|
| Wide Field | 10X | Φ18 | 24.94mm |
| Plan Field | 16X | Φ11 | 15.58mm |

4.3 Total Magnification

| | | | | |
|---------------------|-----|------|------|-------|
| Eyepiece | 10× | 10× | 10× | 10× |
| Objective | 4× | 10× | 40× | 100× |
| Total Magnification | 40× | 100× | 400× | 1000× |

5. Outfit

| Component Name | Specification | Quantity | Standard Outfit |
|----------------|--------------------------------|----------|-----------------|
| Main body | Main Standard | 1 | Optional |
| | Double Layers Mechanical Stage | 1 | Optional |
| | Condenser Holder | 1 | Optional |
| Viewing Head | Digital binocular head | 1 | Optional |
| Condenser | Abbe Condenser, NA=1.25 | 1 | Optional |
| Nosepiece | Quadruple | 1 | Optional |

| | | | |
|--------------|---|-------|----------|
| Illumination | LED illumination | 1 | Optional |
| | 6V20W Halogen Lamp | 1 | Optional |
| | Spare lamp (6V20W Halogen lamp) | 2 | Optional |
| | Spare fuse(2A) | 2 | Optional |
| Eyepieces | 10×Plan Eyepieces | 2 | Optional |
| Objectives | Achromatic objective 4× | 1 | Optional |
| | Achromatic objective 10× | 1 | Optional |
| | Achromatic objective 40× | 1 | Optional |
| | Achromatic objective 100×(oil, spring) | 1 | Optional |
| Condenser | Bright Field Condenser with Adjustable Iris Diaphragm | 1 | Optional |
| Filter | Blue, Green | 1 ea. | Optional |

6. Troubleshooting Guide

1. Optical system

| TROUBLE | CAUSE | SOLUTION |
|--|---|---|
| The edge of the field of view is dark or the brightness is not uniform | The nosepiece is not in the located position (objective and light path not coaxial) | Locate the nosepiece properly where it clicks |
| | The image of filament is not centered | Center the filament |
| | A lens (the objective, condenser, eyepiece or collector) is dirty. | Clean it thoroughly |
| Find dust and stain in the field of view | There are stains on the lens (including condenser, objective, eyepiece and collector) | Clean it up |
| | There are stains on the specimen | Clean it up |
| | The position of the condenser is too low | Loosen the condenser's locking bolt, adjust the condenser to the right position |
| The image is defocused (low resolution \ contrast) | There is no cover slip on the specimen | Add coverslip |
| | The cover slip is too thick or too thin | Use the standard coverslip (0.17mm) |
| | The specimen is placed inversely | Reversal it back |
| | There was oil on the dry objective(easily happened in 40X objective) | Clean it up |

| | | |
|---|---|---|
| The image is defocused (low resolution \ contrast) | There are stains on the lens (including condenser, objective, eyepiece and collector) | Clean it up |
| | didn't use oil for the oil objective | Use immerse oil |
| | There was bleb in the oil | Eliminate the bleb |
| | Use a unsuitable oil | Change to the specified one |
| | The size of the aperture diaphragm is too big | Minify it |
| | There are stains on the incident lens of the binocular tube | Clean it up |
| | The size of the aperture diaphragm is too small | Open it up |
| | The position of the condenser is too low | Adjust the position |
| One side of the image is dark | The condenser is not in the center of the field of view\the condenser inclines | Install the condenser again and adjust the center carefully by centering the bolt |
| | The nosepiece is not in the right position | Turning it until it reach the "clicked" position |
| | The specimen is floating | Fix it |
| The image shift during focusing | The specimen slips on the stage | Fix it |
| | The nosepiece is not in the right position | Turn it to the " clicked "position |
| The image is a little yellow | Not use the blue color filter | Use the blue filter |
| The brightness is not enough | The size of the aperture diaphragm is too small | Adjust again |
| | The position of the condenser is too low | Adjust the position |
| | There are stains on the lens (including condenser, objective, eyepiece and collector) | Clean it up |

2. Mechanical system

| TROUBLE | CAUSE | SOLUTION |
|--|--|--|
| The image can not focus when using high magnification objective | The specimen is placed inversely The coverslip is too thick | Turn inversely Use the standard coverslip (0.17 mm) |
| The objective touch the specimen when changed from low magnification to the higher magnification | The specimen is placed inversely The coverslip is too thick | Turn inversely Use the standard coverslip(0.17 mm) |
| The specimen is not easy to move | The specimen holder is not fixed | Fix it |

| | | |
|---------------------------------------|--|--------------------------------|
| The binocular image is not coincident | The interpupillary distance is not correct | Adjust it |
| Eyes are too tired | No diopter adjustment | Adjust the diopter correctly |
| | The brightness is not suitable | Adjust the voltage of the lamp |

3. Electrical system

| TROUBLE | CAUSE | SOLUTION |
|--|---|--|
| The lamp can't light when the switch is turned on | No power | Check the connection of the power cord |
| | The bulb is not inserted | Insert it correctly |
| | The bulb burns out | Replace it |
| The lamp burns out suddenly | Use a substandard lamp The voltage is too high | Use the specified lamp to replace, if the problem is not solved, contact with the service department |
| The brightness is not enough | Use a substandard lamp The voltage is too low | Use the specified lamp increase the voltage |
| The bulb flickers or the brightness is vertiginous | The bulb is going to burn out | Replace it |
| | The bulb is not entirely inserted into the holder | Check and insert it again |